



### Installation and Set-up

**1** Be sure all parts are included; clamp, arm, locking screw, focus microscope holder.

**2** Put Clamp on the side of the table and tighten the clamp. The width of the clamp opening can be adjusted by moving the two bolts to the lower position.



**3** Insert Screw inside the hole of the clamp.



**4** Put Arm on Screw, and tighten the screw from the top, and cover the hole with a cap.



**5** An adjustable screw is on the rear side of the clamp. Tighten the screw to tighten the arm.



## Installation and Set-up Cont.

- Put focusing rack in the front of the Arm. Slide down the holder and tighten the two top screws, then slide up the holder and tighten the two bottom screws.



- The adjustable screw at the end of the Arm is for adjusting the degree of tightness depending on the weight to be put on the front part of the scope holder.



**Attention:** When lower level position is required for the arm, push the arm down to the position first, and then pull up the focusing rack while pushing down the end of the arm to keep the focus holder at level position.

